

Reveal[®] 3-D for Gliadin R5 Procedure

Note: Please read kit instructions completely before performing test.

Questions? Call +44 (0) 1292 525 610.

Extraction: Rinses




1. Add 0.25 mL of sample into an extraction buffer bottle.



2. Secure cap and shake for 1 minute.

Extraction: Swabs

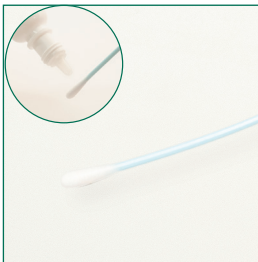

1. Gather the sample. *For dry surfaces moisten with swab wetting solution. Do not premoisten for wet surfaces.*



2. Insert the swab to the extraction buffer bottle and break off into the bottle.



3. Secure cap and shake for 1 minute.

Sample Testing



1. Remove the lid and fill with liquid from the tube.




2. Dip the Reveal 3-D device into the liquid in the lid; ensure that the cavity is saturated with the liquid.



3. Leave the cavity saturated until the liquid is seen running into the test window.



4. Place the device on a flat surface and allow the test to develop for 5 minutes.

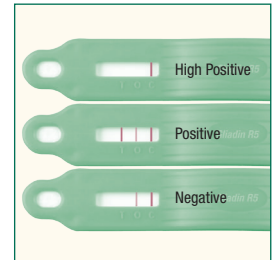

5. Interpret the results.

#8519

Reveal 3-D for Gliadin R5

Rinse: 5–10 ppm
Swab: 2 µg/100 cm²
Testing time: 5 minutes

Materials Provided
20 test devices
20 bottles of extraction buffer
20 sterile swabs

To place an order, or for more information, contact your local Neogen Europe representative, or visit our web site at www.neogeneurope.com

Neogen Europe Ltd. • The Dairy School • Auchincruive, Ayr, Scotland • KA6 5HW
Tel: +44 (0)1292 525610 Fax: +44 (0)1292 525602 • info_uk@neogeneurope.com